

# DIATOME

## Diamond Knives

### Cleaning procedure for our cryo diamond knives


Dear customer,

This procedure serves for the cleaning of our **cryo immuno** and **cryo 25° knives**.

We are at your disposal for any further assistance you might require.



1 Remove the knife stage from the cryo chamber, leave the knives in the stage, wash under tap water for warming up.

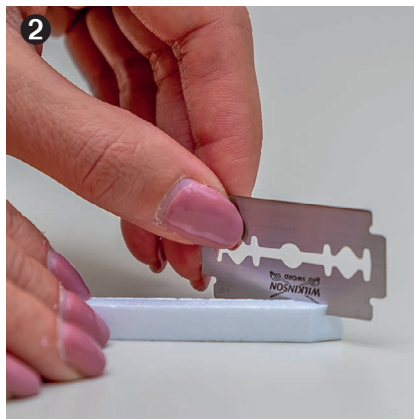

2 Bevel a polystyrene stick with a perfectly clean razor blade.



3 Clean the knife with 15% ethanol, keep wet.



4 Rinse with 100% ethanol.



5 Dry with a dustblower or canned air.

#### DIATOME Ltd

Helmstrasse 1 · CH 2560 Nidau / Switzerland

Phone +41 (0)32 332 91 13

diatome@diatome.ch · www.diatome.ch